

The

# PUREGUARD DISPOSABLE MOPPING SYSTEM



ZERO CAMPAIGN				
INTRODUCTION				
ROTATIONAL CLEANING AND DISINFECTION				
	ANNEX 1 REQUIREMENTS	03		
	DELTA NEUTRAL DETERGENT	04		
PUREGUARD MOPPING SYSTEM				
	MOPS	06		
	HARDWARE	07		
	HOW TO ASSEMBLE THE CLEANROOM TROLLEY	08		
	HOW TO SET UP THE CLEANROOM TROLLEY	10		
	HOW TO ATTACH THE MOP	11		
	HOW TO MOP A CLEANROOM	12		

#### **PUREGUARD BUREGUARD BUREG**

# ZERO

- LIQUID HANDLING
  LIQUID WASTE
  WAITING TIME
  HANDLING OF THE
- MOP SURFACE





### PUREGUARD DISPOSABLE MOPPING SYSTEM

The Micronclean PureGuard disposable mopping system is designed for cleanroom use up to ISO 14644 Class 4, GMP Grade A or B. The system has been designed for ease of use and eliminates the need for

- mixing of chemicals within the cleanroom,
- waiting time within the cleanroom,
- disposal of waste fluids from within the cleanroom and
- the need to disinfect buckets and wringers, which are difficult to clean and sterilise.

In addition, the two-pocket design of the mop enables a hands free and secure connection to the mop frame without the need to handle the mop surface.

Our patented mop system is available in two forms these being:

- Individual mops which are ideally suited to cleanroom floors of less than 150m<sup>2</sup> and
- Multi packed mops which are targeted at cleanroom floors of more than 150m<sup>2</sup>.

The Multi-packed mop option is designed around a mop trolley which features a patented mop pouch dispenser and provides the carrying capacity for both clean and dirty mops to cover the larger floor area.

The heart of the PureGuard system is the pre-saturated microfibre mops which are dosed to give up to 15m<sup>2</sup> coverage. The microfibre mops incorporate extremely low linting 'endless fibre' technology alongside batch manufacturing information to give absolute traceability.

The Micronclean system of pre-saturated mops is fully GMP compliant in that it supports the routine cleaning of the floor using Delta neutral detergent before the application of a rotational disinfection schedule incorporating Alpha and Beta disinfectants.

The Micronclean PureGuard microfibre mops are also available in a dry sterile condition.



# ROTATIONAL CLEANING AND DISINFECTION



## INTRODUCTION

The upcoming revision to EudraLex vol 4 covers Good Manufacturing Practice (GMP) guidelines and Annex 1 is specific to the manufacture of medicinal products. This update comes several years after since its last revision and provides direction to the industry, reflecting best practice.

For cleanroom cleaning and disinfection regimes the key part to note is that the draft guidelines clearly separate the cleaning step from the subsequent disinfection.

<b>☆</b> ☆ ☆ ☆ ☆	☆☆☆☆☆     EUROPEAN COMMISSION ENTERPRISE AND INDUSTRY DIRECTORATE-GENERAL       ☆☆☆☆     Consumer goods Pharmaceuticals
	EudraLex The Rules Governing Medicinal Products in the European Union
	Volume 4
	EU Guidelines to
	Good Manufacturing Practice
	Medicinal Products for Human and Veterinary Use
	Annex 1
	Manufacture of Sterile Medicinal Products
	4 Premises
	4 TTemises
	Disinfection
c	<b>4.36</b> The disinfection of cleanrooms is particularly important. They should be cleaned and disinfected thoroughly in accordance with a written programme. For disinfection to be effective, prior cleaning to remove surface contamination should be performed. More than one type of disinfecting agent should be employed to ensure that where they have different modes of action and their combined usage is effective against all bacteria and fungi.



# **ROTATIONAL CLEANING AND DISINFECTION**

# DELTA NEUTRAL DETERGENT

Micronclean has recently brought to market a brand-new product Delta, which is a neutral detergent. Our Delta product has been developed to be fully compatible with our Alpha and Beta cleanroom disinfection regimes by having the following key features:



# Low foaming formulation

this reduces the risk of streaks and staining when in use which could affect the subsequent disinfectant effectiveness.

#### Neutral (pH 6.5-7.5)

to ensure that it is non-corrosive and will not damage hard surfaces.



### Supplied Pre-mixed and Dosed

to ensure a homogenous distribution of the active ingredient.



### Manufactured with WFI

to ensure maximum cleanroom compatibility.

### Gamma Irradiated to give a sterile product with a two-year shelf life.


# HOW TO USE

With our PureGuard Disposable mopping system, we would recommend the following regime to ensure the maximum benefit from our specially formulated detergent and disinfectants.

Week 1 | Clean the floor by using Delta Mops followed by disinfection using Alpha Mops.

Weeks 2 to 4 | As in Week 1 clean the floor by using Delta Mops followed by disinfection using Alpha Mops.

Week 5 | As in the previous week clean the floor using Delta Mops but then use a Beta Mop as the rotational disinfection step to prevent the potential buildup of resistance within your environment.

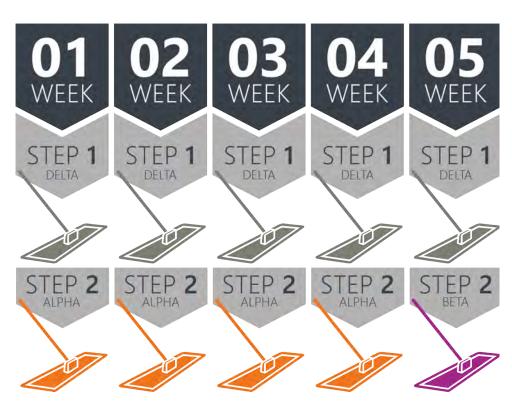



# PUREGUARD MOPPING SYSTEM



# PUREGUARD MOPPING SYSTEM MOPS



### PUREGUARD 1 AND 2 **IPA AND IMS MICROFIBRE MOP** PUREGUARD 3, 4, 5 AND 6 **ALPHA, BETA, DELTA AND DRY MICROFIBRE MOP**

#### FEATURES

Dosed to achieve a uniform coverage of 15m<sup>2</sup>. Proven 2 year shelf life.

Sterile Code	Family	Description	Size	Pack Qty	Case Qty
MSMI01016-A/I	PureGuard 1	IPA Multi Packed Disposable Microfibre Mop Sterile	440 x 140 mm	10	5 x 10
MSMI04016-A/I	PureGuard 1	IPA Individually Packed Disposable Microfibre Mop Sterile	440 x 140 mm	1	40 x 1
MSME01016-A/I	PureGuard 2	IMS Multi Packed Disposable Microfibre Mop Sterile	440 x 140 mm	10	5 x 10
MSME04016-A/I	PureGuard 2	IMS Individually Packed Disposable Microfibre Mop Sterile	440 x 140 mm	1	40 x 1
MSMA01016-A/I	PureGuard 4	Alpha Multi Packed Disposable Microfibre Mop Sterile	440 x 140 mm	10	5 x1 0
MSMA04016-A/I	PureGuard 4	Alpha Individually Packed Disposable Microfibre Mop Sterile	440 x 140 mm	1	40 x 1
MSMB01016-A/I	PureGuard 3	Beta Multi Packed Disposable Microfibre Mop Sterile	440 x 140 mm	10	5 x 10
MSMB04016-A/I	PureGuard 3	Beta Individually Packed Disposable Microfibre Mop Sterile	440 x 140 mm	1	40 x 1
MSMN01016	PureGuard 5	Delta Multi Packed Disposable Microfibre Mop Sterile	440 x 140 mm	10	5 x 10
MSMN04016	PureGuard 5	Delta Individually Packed Disposable Microfibre Mop Sterile	440 x 140 mm	1	40 x 1
MSMD01016	PureGuard 6	Dry Multi Packed Disposable Microfibre Mop Sterile	440 x 140 mm	10	5 x 10
MSMD04016	PureGuard 6	Dry Individually Packed Disposable Microfibre Mop Sterile	440 x 140 mm	1	40 x 1

06 | PUREGUARD DISPOSABLE MOPPING SYSTEM

PUREGUARD MOPPING SYSTEM HARDWARE



## PUREGUARD 7 AND 8 MOP HEADS, TROLLEY AND OTHER ACCESSORIES

Sterile Code	Family	Description	Size	Pack Qty	Case Qty
MTSS01001	PureGuard 7	Mop Trolley	One Size	EA	EA
MFSS01001	PureGuard 7	Lightweight Locking Mop Frame	One Size	EA	EA
MFSS79226	PureGuard 7	Stainless Steel Shrouded Mop Frame (with wing nut)	One Size	EA	EA
MHSS02001	PureGuard 7	PureGuard Stainless Steel Fixed Length Mop Handle (142cm)	One Size	EA	EA
MHSS02002	PureGuard 7	PureGuard Aluminium Adjustable Mop Handle Graphite Grey	One Size	EA	EA
BSP5C1010	PureGuard 8	Polyethylene Bag Sterile	One Size	10	10 x 10

Note: All PureGuard Hardware is fully fully autoclavable.


PUREGUARD DISPOSABLE MOPPING SYSTEM | 07




# **PUREGUARD MOPPING SYSTEM** HOW TO ASSEMBLE A CLEANROOM MOP TROLLEY

Contents

1

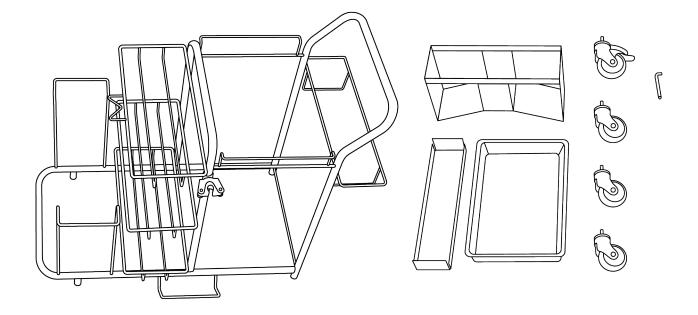




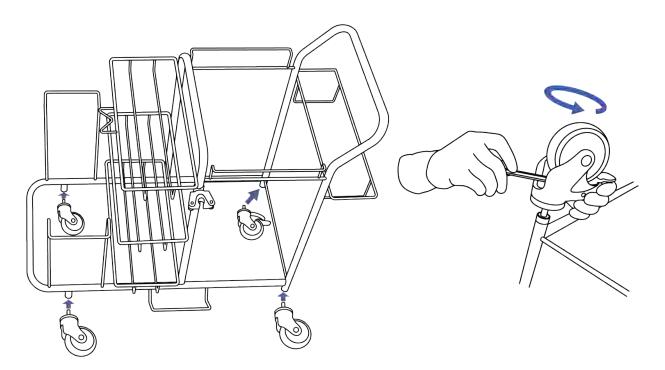

# **PUREGUARD MOPPING SYSTEM** HOW TO ASSEMBLE A CLEANROOM MOP TROLLEY




**3** Positioning the pieces

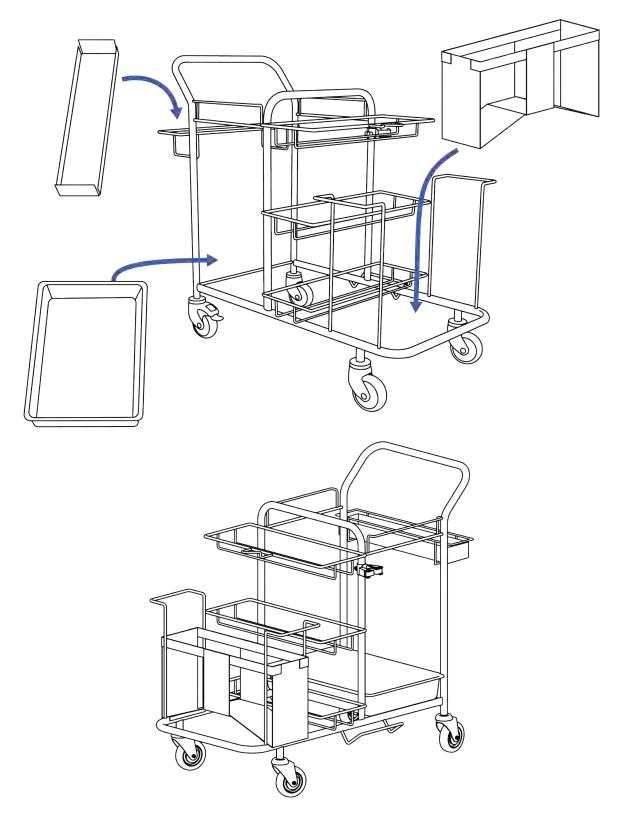



# **PUREGUARD MOPPING SYSTEM** HOW TO SET UP A CLEANROOM MOP TROLLEY

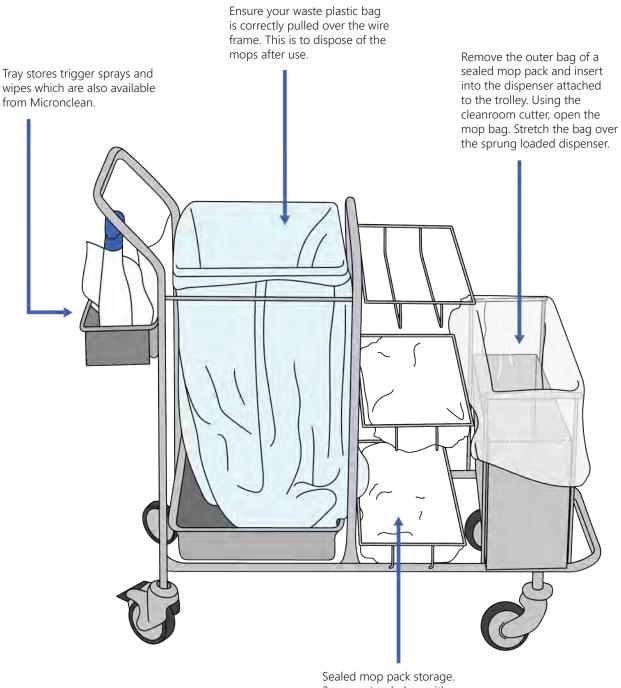


3 open wire shelves with a capacity of 60 mops.

# **PUREGUARD MOPPING SYSTEM** HOW TO USE PUREGUARD MOPPING SYSTEM





Squeeze the metal frame together.

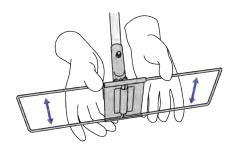



Put the mop frame into the mop dispenser on the front of the trolley.



5

Push the frame into the pockets. Ensure that the mop frame clicks into place.




7

Once the mop is used, squeeze the frame as before over the net bag. Avoid touching the mop head.




2

Push the two sides together.

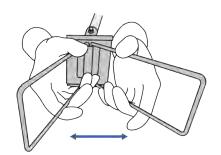



Push down onto the mop head so that the two sides of the frame separate towards the end pockets.

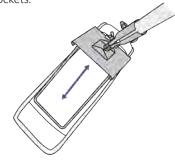



Mop the floor/walls in accordance with aseptic techniques. A single mop cleans up to 15m<sup>2</sup>.



8

Dispose of the mop into the bag. Repeat the process with your next mop. Dispose of the waste bag in accordance with local procedures.







# PUREGUARD MOPPING SYSTEM HOW TO MOP A CLEANROOM

#### The s-curve method

• Mop the two edges either side of your cleanroom area. • To reduce contamination, avoid stepping on any cleaned areas. • Mop across your cleanroom area using an s-shape. Keep lines as neat as possible.

• Overlap paths by 30%.

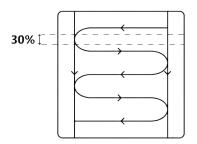

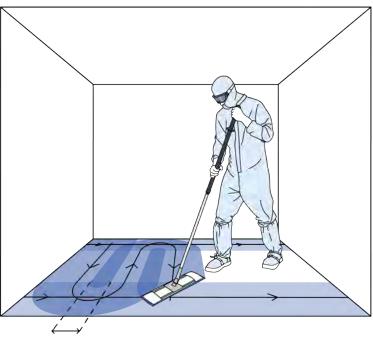

30% overlap

### The lift and lay method


• Place your mop at the edge of your cleanroom area and mop across a straight line.


• Repeat the straight lines, ensuring paths overlap by 30%.

• To reduce contamination, avoid

stepping on any cleaned areas.

• Finally mop along the edge of the room.

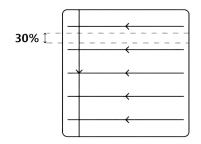

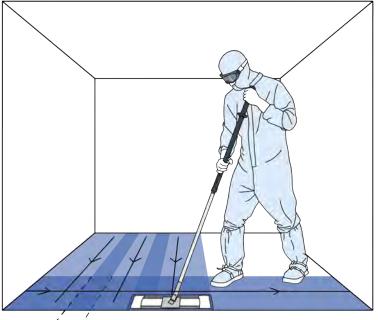

4 30% overlap





Roman Bank / Skegness / Lincolnshire / PE25 1SQ / UK T: +44 (0)1754 767 377 E: customer.services@micronclean.co.uk W: www.micronclean.com