

# INDIVIDUALLY PACKED AUTOCLAVABLE GOGGLES NON-STERILE

Product Code | ONACL0112



VisionGuard **1**

## AUTOCLAVE LOADING PROCEDURE

The Micronclean Autoclavable Goggle has been designed to give outstanding performance in use. To ensure this is achieved in practice Micronclean advise following the autoclave loading method detailed here to reduce the possibility of marks forming on the lenses of the goggles.

### STEP 1 - Sizing of the Autoclave Bag

The correct size of autoclave bag will give enough space between the goggle lens and plastic part of the bag. This minimises the possibility of condensation forming between the goggle lens and the bag.


### STEP 2 - Loading the Goggles into the Autoclave Bag

The autoclavable goggles should always be placed in the selected autoclave compatible bag with the lens upright to prevent any condensation pooling on the lens.



### STEP 3 - Loading the Autoclave

The prepared Autoclave bags should be loaded into the autoclave so that the "paper" is facing down. This again reduces the possibility of condensation pooling on the lens.

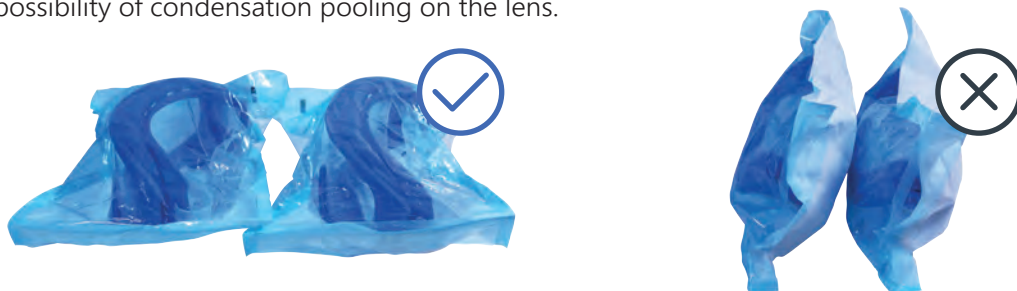

**Please Note:**

1. The Micronclean Autoclavable Goggles are supplied in a Non-Sterile condition within an autoclave bag ready for first processing.
2. The steam used within the autoclave should be 'dry' to avoid problems with drying the goggles during the normal cycle.
3. Do not remove the goggles before the end of the drying cycle.
4. All autoclave procedures should be verified at a local level to ensure they are suitable for purpose.