



# PRODUCT SPECIFICATION

## BETA SESSIONAL POLYCELLULOSE C-FOLDED POUCH WIPES STERILE



### PRODUCT INFORMATION

#### PRODUCT DESCRIPTION

Beta Pre-saturated non-woven Polycellulose sterile wipes.  
Supplied C-folded in a pouch ready immediate use.  
Increased disinfectant and retention from re-sealable pouch.  
Specifically designed for use in Cleanrooms.

#### PRODUCT FEATURES

Non-woven polycellulose wipe with sealed edges.  
Pre-saturated with Alpha to ensure optimum disinfection.  
Supplied in a small pack size to minimise waste.

### TECHNICAL INFORMATION

#### MANUFACTURING

The product is filled and sealed in a Grade B (in use) cleanroom. Packaging is performed in a Grade C (in use) cleanroom.  
The disinfectant is filtered through a 0.2 micron filter at the point of fill.

#### PACKAGING

Supplied double bagged.  
Secondary Packaging: Cardboard carton.

#### STERILISATION DETAILS

Sterilised by Gamma Irradiation (minimum 25kGy).

#### SHELF LIFE

Two years from date of manufacture.

#### STORAGE CONDITIONS

Store in a cool dry place (< 40°C) away from direct sunlight and fluorescent light.

## PRODUCT INFORMATION

### DIMENSIONS

Wipe Size: 23cm x 23cm (9" x 9")

### DISINFECTANT PERFORMANCE

SUSPENSION TESTS							
	Protocol		Organism	Contact Time (min)	Conditions	Log reduction to pass	Log reduction achieved
Suspension Tests Phase 2 Step 1	EN 1276	Bactericidal	Salmonella typhimurium	1 min	Dirty	> 5.0	6.30
			Escherichia coli	1 min	Dirty	> 5.0	6.20
			Enterococcus faecalis	1 min	Dirty	> 5.0	7.30
			Enterococcus hirae	1 min	Dirty	> 5.0	6.60
	EN 1650	Fungicidal/ Yeasticidal	Aspergillus niger	2 mins	Clean	> 4.0	> 4.00
			Candida albicans	2 mins	Clean	> 4.0	> 4.00
	EN 13704	Sporicidal	Clostridium difficile spores	2 mins	Clean	> 3.0	> 3.00
Suspension Tests Phase 2 Step 2	EN 13697	Bactericidal	Pseudomonas aeruginosa	5 mins	Clean	> 4.0	5.70
			Escherichia coli	5 mins	Clean	> 4.0	> 6.56
			Staphylococcus aureus	5 mins	Clean	> 4.0	6.02
			Enterococcus hirae	5 mins	Clean	> 4.0	4.38
		Fungicidal	Aspergillus niger	15 mins	Dirty	> 3.0	4.08

## ORDERING/PACKING INFORMATION

Non-Sterile Code	Sterile Code	Description	Size	Pack Qty	Pack Qty	Case Qty
	WSBT02009	Beta Sessional Polycellulose C-folded Pouch Wipes Sterile	230 x 230mm	15 wipes per Pouch	15	10 x 15

The above product table does not show the regional product codes for clarity. However, to ensure that product labels with the correct languages are ordered and dispatched please contact us.