



PRODUCT SPECIFICATION

IPA POLYCELLULOSE C-FOLDED POUCH WIPE STERILE AND NON-STERILE



PRODUCT INFORMATION

PRODUCT DESCRIPTION

Sterile and Non-sterile nonwoven polycellulose pouch wipe pre-saturated with IPA.

PRODUCT FEATURES

Nonwoven polycellulose wipe.
Supplied C-Folded for ease of use within the cleanroom.
Different packs sizes meet customer requirements.
Pouch is resealable to reduce wastage and improve alcohol retention.

TECHNICAL INFORMATION

MANUFACTURING

The product is filled and sealed in a Grade B (in use) cleanroom.
Packaging is performed in a Grade C (in use) cleanroom.
The disinfectant is filtered through a 0.2 micron filter at the point of fill.

DISINFECTANT SPECIFICATION

Contains: 70 v/v % Propan-2-ol (#67-63-0) and WFI Water for Injection (WFI) grade water.
Unless otherwise stated all Micronclean product labelling for IPA disinfectant products will refer to v/v % as this is the most accurate way of measuring alcohol and water solutions:

	% Tolerance	Notes
IPA 70.0% v/v	+/- 2.0%	@20°C
IPA 62.8% w/w	+/- 2.2%	@20°C

PACKAGING

Primary packaging: Double bagged in polythene packaging.
Secondary Packaging: Cardboard carton.

STERILISATION DETAILS

Gamma Irradiation (minimum 25kGy) sterile variant only.

QUALITY ASSURANCE

Product supplied under BS EN ISO 9001 registered systems.

SHELF LIFE

2 years from date of manufacture.
Use on day of opening.

STORAGE CONDITIONS

Store in cool, well ventilated area.
Keep away from sources of ignition.

MATERIALS & PERFORMANCE

WIPE MATERIAL

55% cellulose & 45% polyester hydroentangled

Performance Data		
Weight	Size	Thickness
68±4g/m ²	228mm x 228mm	0.27mm – 0.34mm

PRODUCT INFORMATION

DISINFECTANT PERFORMANCE

SURFACE TESTS						
Protocol	Organism	ATCC Code	Condition	Log reduction to pass	Log reduction @ 5 mins	Outcome
BS EN 13697:2015	S. aureus	6538	Clean	Log 4.0	> 6.62	Bactericidal
	P. aeruginosa	15442	Clean	Log 4.0	> 6.83	Bactericidal
	E. coli	10536	Clean	Log 4.0	> 6.85	Bactericidal
	E. hirae	10541	Clean	Log 4.0	> 6.66	Bactericidal

SURFACE TESTS						
Protocol	Organism	ATCC Code	Condition	Log reduction to pass	Log reduction @ 15 mins	Outcome
BS EN 13697:2015	C. albicans	10231	Clean	Log 3.0	> 5.26	Yesticidal
	A. brasiliensis	16404	Clean	Log 3.0	3.04	Fungicidal

Please note: Data shown above are just typical values.

ORDERING/PACKING INFORMATION

Non-Sterile Code	Sterile Code	Description	Size	Pack Qty	Case Qty	Case Dimensions	Weight
WNVG04023		IPA Polycellulose C-folded Pouch Wipe Non-sterile	230x230mm	100	8 x 100	310 x 280 x 390mm	8.5 Kg
	WSVG04023	IPA Polycellulose C-folded Pouch Wipes Sterile	230x230mm	100	8 x 100	310 x 280 x 390mm	8.5 Kg
	WSCW92009	IPA Sessional Polycellulose C-folded Pouch Wipe Sterile	230x230mm	15	40 x 15	405 x 405 x 195mm	8.44 Kg

The above product table does not show the regional product codes for clarity. However, to ensure that product labels with the correct languages are ordered and dispatched please contact us.