



PRODUCT SPECIFICATION

IMS SILVER PACK

POLYETHYLENE/POLYPROPYLENE

Z-FOLDED WIPES STERILE



PRODUCT INFORMATION

PRODUCT DESCRIPTION

VeriGuard 9 is pre-saturated pouch wipe designed for use in lower grade cleanrooms as a hard surface wipe. The resealable pouch gives increased alcohol retention.

PRODUCT FEATURES

Non-woven Polyethylene/Polypropylene Z-folded sealed edge wipes. Pre-saturated with IMS 70 /30 DI. Z-Folded within the re-sealable pouch.

MANUFACTURING

Packed and pre-saturated under controlled conditions.

DISINFECTANT SPECIFICATION

Contains: 70 v/v % IMS (Denatured alcohol Ethanol and Propan-2-ol) and De-Ionised Water (DI) grade water.

Unless otherwise stated all Micronclean product labelling for IMS disinfectant products will refer to v/v % as this is the most accurate way of measuring alcohol and water solutions:

	v/v %	CAS No	% Tolerance	Notes
Ethanol (ethyl alcohol)	66.5	64-17-5	+/- 2.0%	@20°C
Propan-2-ol	3.5	67-63-0	+/- 2.0%	@20°C

PACKAGING

Packed in cleanroom compatible pouches containing 200 wipes. Double Bagged.

STERILISATION DETAILS

Sterilised by Gamma Irradiation (minimum 25kGy).

QUALITY ASSURANCE

Product supplied under BS EN ISO 9001.

SHELF LIFE

2 years from date of manufacture.

STORAGE CONDITIONS

Store in a cool, dry, well ventilated area and protect from direct sunlight heat and naked flames.

MATERIALS & PERFORMANCE

Performance Data

Weight

23g/m² +/-5%

Size

220mm x 200mm (9" x 8")

PRODUCT INFORMATION

DISINFECTANT PERFORMANCE

SURFACE TESTS

Protocol	Organism	Code	Condition	Log reduction to pass	Log reduction @ 5 mins	Outcome
BS EN 13697:2015	P. aeruginosa	DSM 939	Clean	Log 4.0	> 6.57	Bactericidal
	E. coli	NCTC 10418	Clean	Log 4.0	> 6.23	Bactericidal
	S. aureus	NCTC 10788	Clean	Log 4.0	> 6.44	Bactericidal
	E. hirae	DSM 3320	Clean	Log 4.0	> 4.41	Bactericidal
Protocol	Organism	Code	Condition	Log reduction to pass	Log reduction @ 15 mins	Outcome
BS EN 13697:2015	C. albicans	DSM 1386	Clean	Log 3.0	> 6.29	Yesticidal

SUSPENSION TESTS

Protocol	Organism	Code	Condition	Log reduction to pass	Log reduction @ 5 mins	Outcome
BS EN 1276:2015	P. aeruginosa	DSM 939	Clean	Log 5.0	> 5.26	Bactericidal
	E. coli	NCTC 10418	Clean	Log 5.0	> 5.54	Bactericidal
	S. aureus	NCTC 10788	Clean	Log 5.0	> 5.17	Bactericidal
	E. hirae	DSM 3320	Clean	Log 5.0	> 5.17	Bactericidal
Protocol	Organism	Code	Condition	Log reduction to pass	Log reduction @ 15 mins	Outcome
BS EN 1650:2015	C. albicans	DSM 1386	Clean	Log 4.0	> 4.26	Yesticidal

Please note: Data shown above are just typical values.

ORDERING/PACKING INFORMATION

Sterile Code	Description	Size	Pack Qty	Case Qty	Case Dimensions	Weight
WSCW97ZZZ	VeriGuard 9 IMS Silver Pack Polyethylene/Polypropylene Z-folded Wipes Sterile	200mm x 220mm (8" x 9")	200	8	530 x 235 x 190mm	7.5kg

The above product table does not show the regional product codes for clarity. However, to ensure that product labels with the correct languages are ordered and dispatched please contact us.